**Tris-NaCl Buffer,
with Albumin
pH 7.40
Set of 4x20 mL**

REF. AR005L

For *in-vitro* use only

PRESENTATION:

Buffer containing Tris (Tris (hydroxymethyl) aminomethane) (6.1 g/L), Sodium Chloride (8.77 g/L), Bovine Serum Albumin (BSA at 10 g/L) and 0.9 g/L Sodium Azide (as preservative).

*CAUTION: Sodium Azide (NaN₃) may react with lead and copper plumbing to form highly explosive metal azides. Flush with large volumes of water when discarding into a sink.

This kit is for research use only and should not be used for patient diagnosis or treatment.

USE:

For *in vitro* use only.

Vials of 20 mL, ready to use.

Dilution buffer for biological tests, recommended mainly for diluting purified proteins and enzymes (Albumin prevents adsorption of purified proteins on recipient walls).

Can be used as buffer recommended by the US and EU Pharmacopoeia for the assay of heparin anti-IIa and anti-Xa activity, verified and validated in combination with BIOPHEN™ ANTI-IIa (2 Stages Heparin Assay, ref. 220005) and BIOPHEN™ ANTI-Xa (2 Stages Heparin Assay, ref. 221005) kits.

Notes:

- For instructions concerning the tests to be performed, refer to the specific associated kit inserts.
- Take care of stabilizing the solution at room temperature (18-25°C) prior to use.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

PH:

7.40

STABILITY:

In its original vial, unopened and kept at 2-8°C, stable until the expiration date printed on the label.

When open, protected from any contamination, this solution is stable for 4 weeks at 2-8°C.

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